

Information	Page 1 of 2
Shipping of MBASS30 for Maintenance & Recalibration	Date: 01.10.2024 Revision: 2 Created: hh

Document: Infoblatt_Verpackung_MBASS30_en_5-3-2-2-11_2_20241001.docx

Scope of Delivery

For maintenance and recalibration, we **only** need the parts listed below – the mentioned sampling heads as available:

- MBASS30
- Charger with power cord
- Sampling head LKS 30
- Sampling head LKS100
- Particle sampling head PS 30 with slide holding plate
- Filter adapter FA 30
- Completed Declaration of Contamination (<https://www.holbach.biz/en/Downloads>)
- Handling case

Notes:

- The charger gets only tested for the provided charging current. We do not perform an electrical safety check.
- We use the provided sampling heads of the customer for the recalibration. The serial numbers of the sampling heads will be specified in the calibration certificate.
- **Prior** to shipping, remove all parts that are not required, such as: cables, adapters, adhesive tape, disinfectants, user manuals, calibration certificates, scissors, tweezers and other tools. Unneeded accessories increase the transport weight and thus the shipping costs. Accessories and tools lying loosely in the handling case may damage the device during transport.

Packaging

Ideally, MBASS30 and the accessories mentioned above are shipped to us (in the handling case) in the original packaging.

If the original packaging is no longer available, please use a sturdy (double wall) cardboard box. Protect the parts sufficiently with bubble wrap or similar dust-free padding material.

Suitable for dust-free padding of the device and the accessories are:

- ✓ Bubble wrap
- ✓ Air cushion pads
- ✓ Foam material

The following padding materials are **not suitable**, because of possible dust:

- ✗ Paper
- ✗ Packing peanuts (also compostable ones)
- ✗ Expanded polystyrene
- ✗ Wood wool

Note: If no dust-free padding material is available, please wrap the parts in plastic bags (e.g. a new trash bag) and seal the bags carefully.

Information	Page 2 of 2
Shipping of MBASS30 for Maintenance & Recalibration	Date: 01.10.2024 Revision: 2 Created: hh

Document: Infoblatt_Verpackung_MBASS30_en_5-3-2-2-11_2_20241001.docx

We will return the device and the components in the same packaging in which we received them from you. We do not assess whether the packaging is appropriate and suitable for the device. We are therefore not liable for transport damage due to unsuitable packaging. We are expressly not liable for damage or loss both when receiving the devices and when returning them.

Notes on shipping in the handling case:

- ✓ Put a sampling head onto the MBASS30 system. This protects the locking springs and closes the sampling air inlet.
- ✓ Pack the upper part and the slide holding plate of the particle sampler PS 30 together in one bag. That way, the holding plate cannot fly around in the handling case.
- ✓ Pad the parts in the handling case on top with dust-free padding material (see above).



Figure 1: Parts for shipping prior to padding

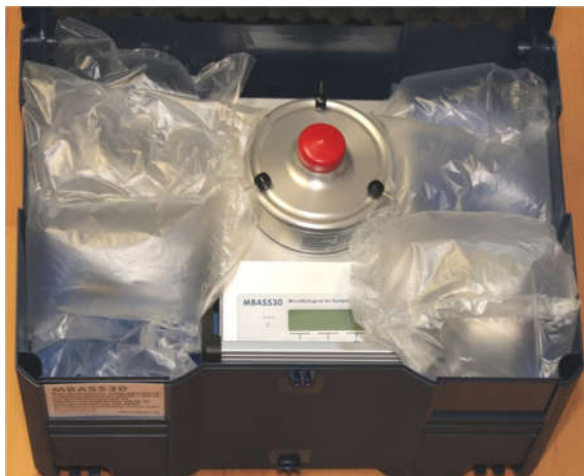


Figure 2: Padded accessories

Please do not do this:



Figure 3: Situation as delivered without padding and with unnecessary accessories



Figure 4: unnecessary damage of the side panel caused by missing padding during transport